

DOWNLOADING FLUXERGY WORKS & SAMPLE TEST IMPORT

Downloading Fluxergy Works

Download link: <https://fluxergy.com/downloads/Fluxergy-Works-CE.exe>

Serial Number: Ask your Fluxergy Rep for a Serial number to activate your software

A couple of things to note here when installing Fluxergy Works:

Enter the serial number provided to you by Fluxergy.

- a. Note: there is a one-time activation process which **requires an internet connection**. After activation internet access is no longer necessary.

After installing software, create an administrator account.

- a. Note: the administrator account created during Fluxergy Works install is the master administrator and cannot be edited by other administrator accounts.
- b. Note: **if the master administrative account password is lost, it cannot be reset**. To mitigate the risk of losing all administrative access, Fluxergy recommends creating a secondary administrator account immediately after install
- c. If all user admin passwords are lost, users must uninstall and reinstall Fluxergy Works after backup of test files and deletion of test storage folders from previous installs of Fluxergy Works

Ask your Fluxergy rep for a download of an .flx file (specific file type for Fluxergy Works) that you can import into Fluxergy Works. This will have an example test to show your users!

Simply select import in Fluxergy Works on the history tab and select the file. You can view the sample test.

History Search Here

History Imported

Test ID	Test Name	Template Name	Device Name	Status	Date
2	VT001	CoVID-19 Sample-to-Answer-RT-...	192.168.10.4	Completed	2020-09-29 17:35:54.662
1	denis	CoVID-19 Sample-to-Answer-RT-...	192.168.0.16	Completed	2021-03-17 06:51:31.763

IMPORT

Dashboard

Template

Run Test

History

Device

Admin

Logout